

PRODUCT DESCRIPTION

This kit is to be used to replace the tank discharge hose on an Encon Gravity-Fed Eyewash.

FIGURE 1

NORMAL OPERATING POSITION VIEW



PIVOT PIN LOCATION

NOTICE:

Modifications to the Encon Gravity-Fed Eyewash tray and closure mechanism require that you identify the style unit you have before proceeding to replace the tank discharge hose. The Gravity-Fed Eyewashes built prior to April 1990 CANNOT be repaired. Only units whose trays are held closed by means of an interlocking male/female black latch mechanism, as shown in **Figure 1**, are repairable. Replacement units can be ordered from your local Encon distributor.

COMPONENT PARTS LIST

Kit includes one (1) 7" Hose and two (2) Hose Clamps.

INSTALLATION INSTRUCTIONS

1. If tray does have a black interlocking male/female latch mechanism, examine the upper tray section.
 - a) If the "U" canal hose path is 7/8" depth (see **Figure 2**), then proceed with hose replacement.
 - b) If "U" canal hose path is 1/2" or less (see **Figure 3**), an upper tray replacement section is required (P/N 01110711).
2. Pull down tray to normal operating position as shown in **Figure 1**. Remove four (4) screws from the upper tray section and remove lower tray section. Remove pivot pins and set aside. Lay tank on its back for ease of hose replacement.
3. Remove and discard old discharge hose by removing hose clamps from barbed fittings.
4. Place two (2) new hose clamps on new 7" hose provided. Push one end of the hose onto barbed fitting until flushed with tank. Position and tighten clamp on barbed fitting with needle nose pliers for water tight connection. Push other end of hose completely onto the barbed fitting of the tee in the tray top and secure with second clamp.
5. Lay upper tray section in the closed position with the new hose stretched in the center "U" canal hose path at the rear of the tray. Replace pivot pins and snap into position with the chamfered end inserted into the tank pivot cavity (see **Figure 1**).
6. Lay lower tray section into position on upper tray section. While squeezing both sections, lower tray assembly to normal operating position, exercising care to prevent the separation of the tray halves, and secure screws/nuts. Inspect the clamp on hose to ensure hose is completely secured on barbed fitting.
7. Refill eyewash station with clean potable water and Hydrosep® water preservative. Follow directions on Hydrosep® water preservative container.

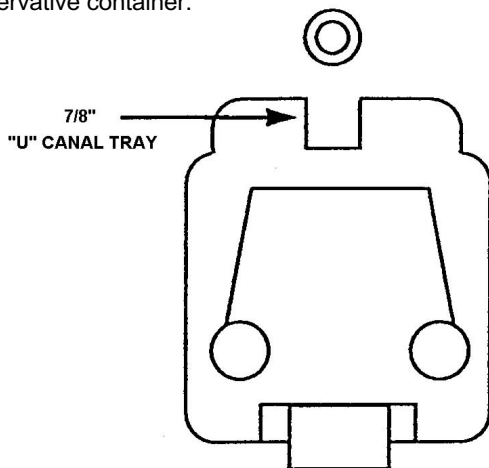


FIGURE 2

NEW TRAY

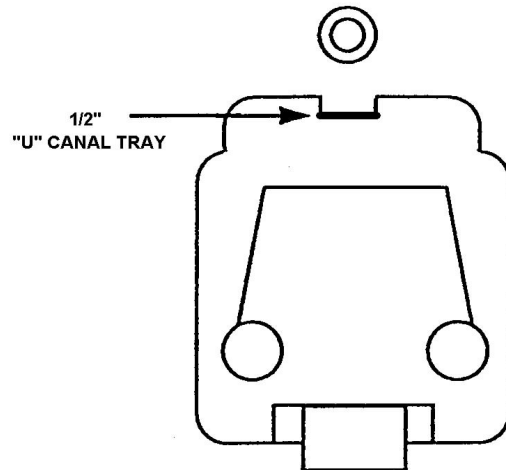


FIGURE 3

OLD TRAY

PROPOSITION 65

WARNING: This product can expose you to BPA, which is known to the State of California to cause birth defects or other reproductive harm. For more information go to www.P65Warnings.ca.gov.

WARRANTY STATEMENT

ENCON HEREBY DISCLAIMS ALL WARRANTIES EXPRESSED OR IMPLIED INCLUDING BUT NOT LIMITED TO WARRANTIES OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, AND NON-INFRINGEMENT OF THIRD-PARTY RIGHTS, EXCEPT AS HEREINAFTER PROVIDED.

Encon Safety Products warrants that for one year from the date of purchase of any Encon products, the product will be free of defects in materials and workmanship if properly used and cared for or cleaned under normal conditions in accordance with Encon's use and care instructions and properly installed, if applicable, in accordance with Encon's installation instructions. With respect to the product, Encon's only obligation and purchaser's exclusive remedy under this warranty is to repair or replace such product; provided that:

1. Encon is notified of the defect within one year of shipment, and
2. the product is determined by Encon to be defective.

Encon requires proof of original ownership as proof of warranty coverage, and Encon must receive any claim under this Limited Warranty within one year of purchase of the product.

NOTWITHSTANDING ANYTHING TO THE CONTRARY CONTAINED HEREIN, ENCON SHALL NOT BE LIABLE FOR LOSS, DAMAGE, OR EXPENSE ARISING DIRECTLY OR INDIRECTLY AS A CONSEQUENCE OF USE OF THE EQUIPMENT WITH OTHER PRODUCTS OR FROM ANY OTHER CAUSE, INCLUDING ANY CONSEQUENTIAL, INCIDENTAL, SPECIAL OR EXEMPLARY DAMAGES, EXCEPT FOR ENCON'S OBLIGATION TO REPAIR OR REPLACE DEFECTIVE PRODUCTS AS EXPRESSLY PROVIDED IN THIS STATEMENT.

Replacement parts purchased from Encon are warranted for one year following the shipment of such replacement part, or until the expiration of the warranty period for the product, whichever is less. No warranty is given in connection with products that are altered without Encon's expressed written consent. The same warranty limitations and the obligations of Encon as set out herein above shall apply to replacement parts.

Encon's total liability arising out of this warranty (including, but not limited to, warranty claims) regardless of forum and regardless of whether such action or claim is based on tort, contract or otherwise will not exceed the total purchase price of the product.



6825 W. Sam Houston Pkwy. N
Houston, TX 77041
1-800-AT-ENCON / 1-800-283-6266
customerservice@enconsafety.com
<http://www.enconsafety.com>